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042-719-4999 / 090-6171-7744

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PRODUCT MANUAL

FOR THE

ANION SELF-REGENERATING SUPPRESSOR 300
(ASRS[®] 300 (4-mm), P/N 064554)
(ASRS[®] 300 (2-mm), P/N 064555)

CATION SELF-REGENERATING SUPPRESSOR 300
(CSRS[®] 300 (4-mm), P/N 064556)
(CSRS[®] 300 (2-mm), P/N 064557)

©Dionex Corporation, 2007
Document No. 031956
Revision 05
August 2007

1. INTRODUCTION

1.1. Self-Regenerating Suppressor

The Self-Regenerating Suppressor (SRS® 300) replaces the ULTRA II suppressor product line. The SRS 300 is redesigned with exterior hardware changes that allow the suppressor to be more pressure tolerant than previous generation suppressor devices. The SRS 300 continues to use cleaned ion exchange components (gasketed screens and membranes) as the ULTRA II suppressor devices. The 2mm ASRS 300 replaces the ASRS MS suppressor product and should be used for anion Mass Spectrometry (MS) Applications.

The Self-Regenerating Suppressor 300 (SRS 300) is available in two versions: the Anion Self-Regenerating Suppressor 300 (ASRS® 300) or Cation Self-Regenerating Suppressor 300 (CSRS® 300). The SRS 300 system consists of a Self-Regenerating Suppressor, the SRS Control, the back pressure coils, and the Gas Separator Waste Tube. This high performance, low maintenance AutoSuppression® system provides a reliable solution for Ion Chromatography.

Additionally, the SRS 300 offers high capacity suppression while adding minimal dead volume to the analytical system. The ASRS 300 provides continuous suppression of traditional eluents, and more concentrated eluents up to 200 mM NaOH. The CSRS 300 offers continuous suppression of concentrated eluents up to 100 mN H₂SO₄ or MSA. This high capacity significantly expands the capabilities and simplifies the operation of Ion Chromatography.

The SRS 300 is available in both 2-mm and 4-mm formats for use with 2, 3, 4, or 5-mm Ion Chromatography columns and systems. The 2-mm SRS 300 is specially designed with reduced internal volume to ensure optimum performance with 2-mm columns and systems. In fact, the suppression capacity of the 2-mm version is greater than the 4-mm version due to the reduced eluent flow rate and proportionally greater surface area.

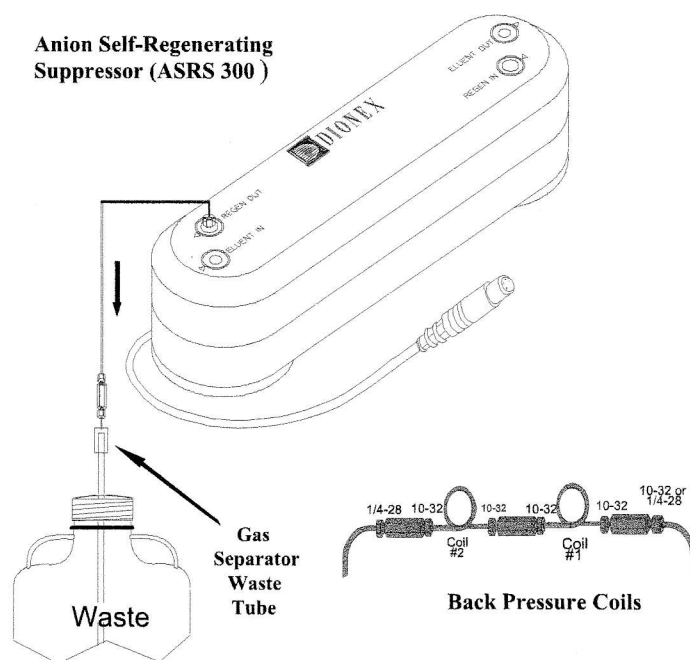
Figure 1
The Self-Regenerating Suppressor

Assistance is available for any problem during the shipment or operation of Dionex products through the Dionex North America Technical Call Center at 1-800-DIONEX-0 (1-800-346-6390) or through any of the Worldwide Dionex Offices on the Reference Library CD-ROM.