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042-719-4999 / 090-6171-7744

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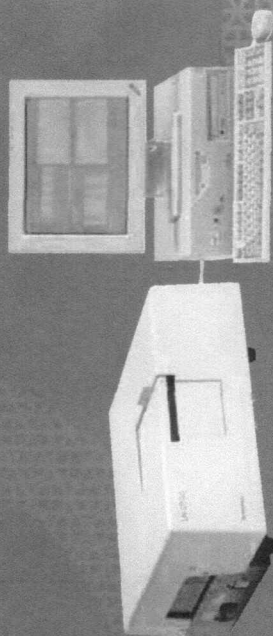
Ultra low stray light in the world's best
double monochromator instrument
UV-2550
High performance
single monochromator instrument
UV-2450

UV-2550 UV-2450

Shimadzu UV-VIS Spectrophotometers

UV-2550 Ultra low stray light in the world's best double monochromator instrument
UV-2450 High performance single monochromator instrument

Shimadzu's highly reputed spectrophotometers combined with the most up-to-date, superb software achieve high throughput and increased productivity. With the newly developed 8/16 cell positioner system, up to 16 specimens can be analyzed automatically.

Low stray light

The UV-2550, with its DDM (double biased grating double monochromator) design, provides a high energy throughput optical system with ultra low stray light. The UV-2450, the single monochromator version of the UV-2550, achieves 0.015% stray light even with its simpler optical system.

High Performance

Both instruments offer outstanding cost performance in their ability to provide UV-VIS spectral analysis of wide ranging samples, including organic and inorganic compounds, biological samples like DNA and enzymes, as well as special measurements on optical materials.

Windows Compatible

Operating under the Windows XP/2000 environment, UVProbe software is provided as standard. UVProbe incorporates the latest multitasking and report generation functionality, including network support, enabling measurement assistance among colleagues and effective presentation of previous data.

Analysis of multiple trace samples

The 8/16 microtitercell permits selection among 4 possible combinations with sample capacities of 25µL and 100µL, and either the normal type or the constant-temperature controlled type cell holder may be selected.

NOTE! Both the cell holder and the multi-cell are optional accessories.

(100 µL multi-cell)

Contents

P 04 - Features

P 12 - Specifications

P 13 - Option

P 10 - Applications

6383

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©CISTEC